

CARSON®

MicroBrite™ Plus

60-120x LED Lighted Pocket Microscope



MM-300

Congratulations on selecting your new MicroBrite™ Plus Microscope! In order to achieve optimum performance, please follow instructions for proper use and care.

PRODUCT SPECIFICATION


Magnification:	60x -120x	Weight:	0.13 lbs (excluding battery)	Battery:	1 AA (not included)
Exit Pupil Diameter:	10mm - 14mm	Dimensions:	2.3" x 1.2 x" 3.7"		
Optical zoom:	2x	Illumination:	LED		

INSTALLING THE BATTERY

1. Slide open battery door.



2. Insert 1 AA Battery with the correct polarity and replace battery door.

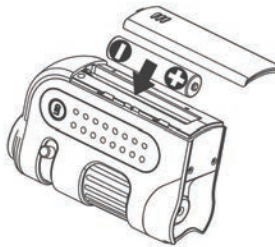

WARNING: Use only alkaline battery. Keep away from fire or excessive heat. Inspect battery for corrosion or leakage. Insert in direction indicated. Rechargeable battery not recommended.

USING THE MICROSCOPE

1. Place viewing object directly below Objective Lens.



2. Press LED switch to turn light On and Off.



3. Adjust Zooming Dial to desired magnification.



4. Rotate Focusing Ring until image is clear and sharp.



WARNING

- To avoid permanent eye damage:
 - Do not look at the sun with the microscope.
 - Do not place the lighted microscope directly upon human eye.
- Remove battery before storing microscope for extended periods.
- The device may get warm after extended use. This is normal and should not be considered a defect.

Customer Service:

If you experience any difficulties, please contact us and we will be happy to help you.



✉ info@carson.com



✉ uksupport@carson.com



✉ eusupport@carson.com

For warranty information, visit www.carson.com/warranty

www.carson.com

Carson Optical 2070 5th Avenue, Ronkonkoma, NY 11779 | Tel: 631-963-5000